


## Viewing chromatograms in the trend data window

It is possible to view the chromatogram that is associated with a particular trend data point.

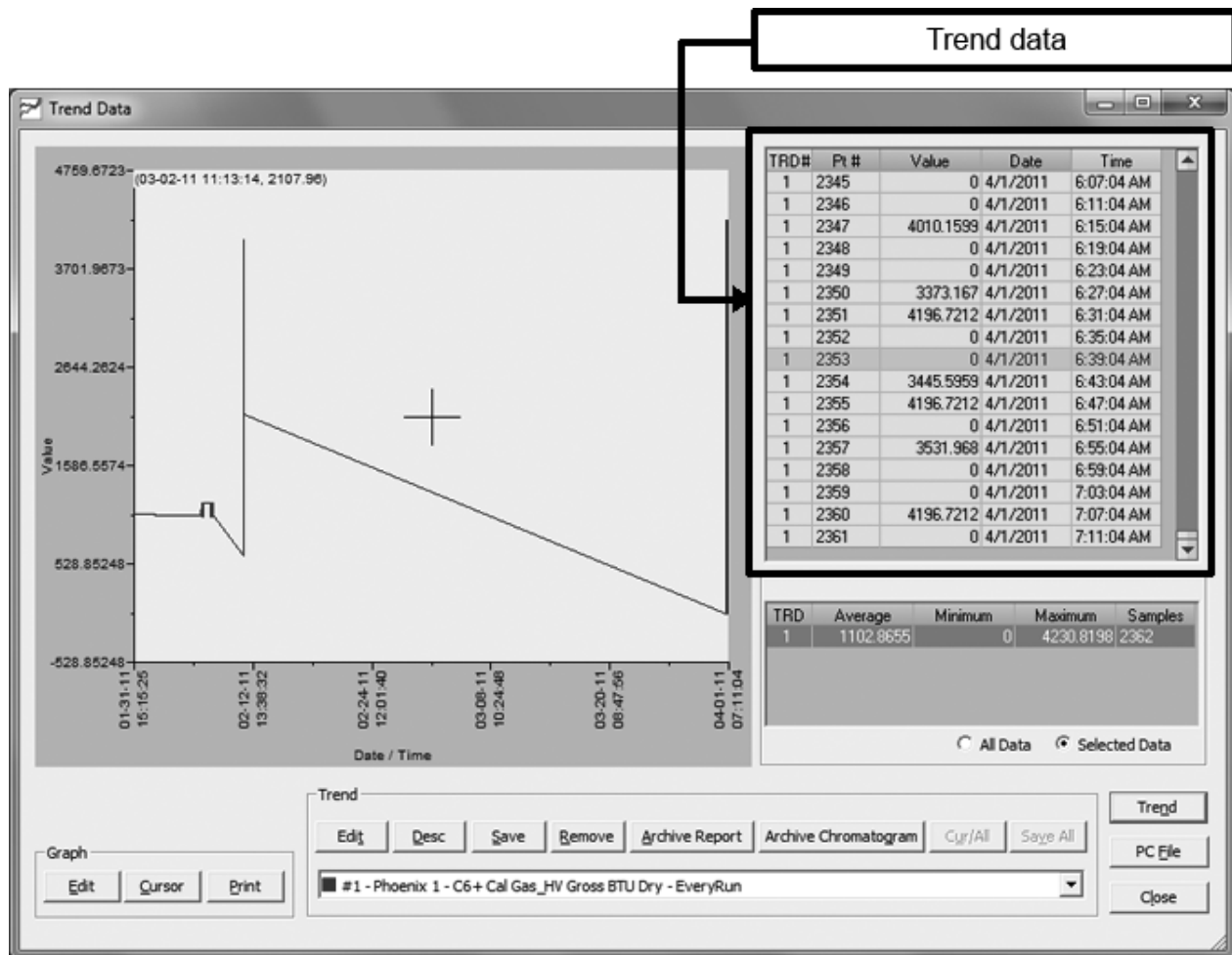

**Note:** This feature is only available for **live** trends.

**Note:** This feature is limited to trends whose data is approximately five days old or less. This is because the GC typically stores a chromatogram for approximately five days before it is deleted to make room for newer chromatograms.

To view an associated chromatogram from the *Trend Data* window, do the following:

1. Select the trend point (Pt. #) from the trend data table.

**Note:** You can use the CTRL or SHIFT key to select more than one trend point.

2. Click **Archive Chromatogram**. The *Chromatogram Viewer* opens with the appropriate chromatogram displayed.

**Note:** If you selected more than one chromatogram, they will all display on the same window, each in a different color.

